

How to find a serial number

- Option 1:

Have a look at the rear side of the instrument. The serial number starts with the letters “MVE”




- Option 2:

Open the settings page of the Phenom user interface.

For legacy user interface (all instruments except XL G2, Pharos G2, Pure G6, Pro G6 and ProX G6)

- Activate the “advanced mode” with the PW “expert”
- Go to “Advanced”, page “Network settings”
- The Instrument ID can be found on the left bottom area

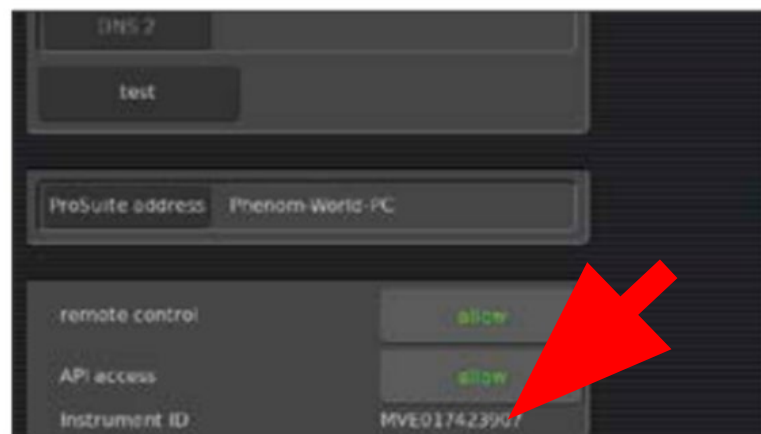

For new user interface (only instruments Phenom XL G2, Pharos G2, Pure G6, Pro G6 and ProX G6)

- Goto settings, page “Status”
- The serial number starts with the letters “MVE”